

Serial Dilutions


Serial dilutions are often performed when titrating antibodies or when generating amplified dilutions of an analyte. A serial dilution is a step-wise and geometric series of dilutions which starts with a small amount of starting material and amplifies the dilution factor serially by using diluted material as a source for subsequent dilutions. Serial two-fold and ten-fold dilutions are commonly used to titer antibodies or prepare diluted analytes (for a standard curve for example). Serial dilutions are also commonly used to avoid having to pipette very small volumes (1-10 μl) to make a dilution of a solution. Serial dilutions can be performed in tubes (microfuge or 12 X 75 mm) or in the wells of plates (depending on the volumes you require). It is important to know what volume of the dilution you will need and the range of concentrations that you desire.

In the example below we are performing two-fold serial dilutions of a $\frac{1.0 \text{ mg}}{\text{ml}}$ stock of analyte for use as standards. We would like the concentration of the standards to be in the range of $\frac{1000 \text{ ng}}{\text{ml}}$ to about $\frac{10 \text{ ng}}{\text{ml}}$. The original or first dilution will dilute the stock for the highest primary concentration that we need, $\frac{1000 \text{ ng}}{\text{ml}}$. This will be done by performing an initial 1: 1000 dilution because:

$$\frac{C1}{C2} = \text{dilution factor}$$

$$\frac{1,000,000 \text{ ng/ml}}{1,000 \text{ ng/ml}} = 1000 \text{ dilution factor}$$

To prepare the initial 1:1,000 dilution we can take 5 μl of the stock solution and add 4,995 μl of diluent (buffer or water). This would result in a solution of analyte with a concentration of $\frac{1000 \text{ ng}}{\text{ml}}$. Now the two-fold dilutions can be prepared from the $\frac{1000 \text{ ng}}{\text{ml}}$ solution. For the first 1:2 dilution we would take 500 μl of the 1:1,000 dilution and add it to 500 μl of diluent in the next tube. This will result in a solution with a concentration of $\frac{500 \text{ ng}}{\text{ml}}$. This solution is then used in a subsequent or serial dilution by taking 500 μl of the previous $\frac{500 \text{ ng}}{\text{ml}}$ solution and adding it to the next tube containing 500 μl of diluent. This dilution will now result in a solution with a concentration of $\frac{250 \text{ ng}}{\text{ml}}$. This process is repeated until a final solution at a concentration near $\frac{10 \text{ ng}}{\text{ml}}$ is made. In the example below, the highest dilution has a concentration of $\frac{7.8 \text{ ng}}{\text{ml}}$.

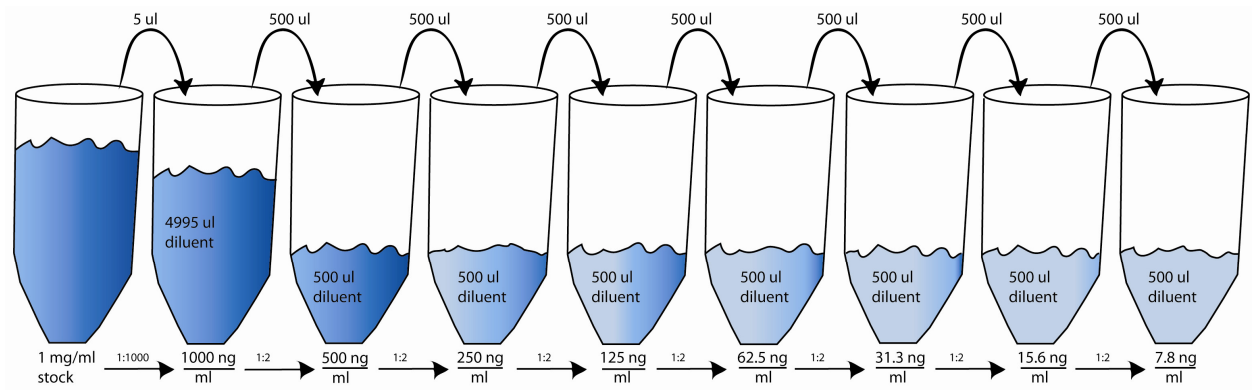

The total dilution factor for any of the resulting solutions can be calculated by multiplying each dilution factor leading up to a particular dilution. For example if we wanted to find the total dilution for the $\frac{62.5ng}{ml}$ solution :

$$DF_{total} = DF1 \times DF2 \times DF3 \times DF4 \times DF5$$

$$DF_{total} = 1000 \times 2 \times 2 \times 2 \times 2$$

$$DF_{total} = 16,000$$